Heating and Cooling from 5 to 85 °C

- Versatile
- High Accuracy
- Even Air Distribution
- Simple Control

KNAUER's **Smartline Column Thermostat** has been specially designed for the use in chemical and biological routine analysis, including research laboratories. Its areas of application range from analytical to preparative HPLC.

A stable column temperature not only improves the reproducibility, but rather considerably improves the resolution of certain HPLC methods as well.

The reliable control of the temperature, which covers a range of 5 - 85 °C has been realized with the use of a microprocessor controlled Peltier element. The measurement of the temperature is carried out with high accuracy using double reference sensor technology.

The powerful interior fan provides for 2-way air circulation and a perfectly even distribution of the thermostatted air in the column compartment, enabling quick ramping of the temperature.

The remote keypad enables access to all of the thermostat's functions and displays target and actual temperature in increments of 0.1 °C.

Depending on the column diameter the **Smartline Column Thermostat** can hold up to four liquid chromatography columns with a maximum length of 350 mm each. The maximum inner diameter of the HPLC column is 20 mm.

In addition to the standard temperature adjustment, it is possible to program temperature gradients with up to 99 steps to optimize an analysis.

The temperature can either be controlled with the supplied control panel or externally with a personal computer. Gradients can be conveniently programmed with our chromatography software ChromGate®.

The **Smartline Column Thermostat** features a power-on self-check, a leak detector with adjustable sensitivity, and a programmable turn-off temperature to protect the column from overheating.

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Technical Data and Ordering Information

Heating and Cooling System

Microprocessor controlled Peltier element for heating and cooling. Fan supported 2-way air circulation

Column Compartment

Interior thermally insulated

Columns analytical and preparative columns up to 20 mm ID and 350 mm length; maximum of 4 columns

Dimensions (interior) 90 x 390 x 47 mm (W x H x D)

Injection Valve optional with an integrated 6-Port/2-channel injection valve (stainless steel or PEEK)

Temperature Control

Temperature Range 5 - 85 °C

Accuracy ± 0.2 °C

Temperature Stability ± 0.1 °C

Heating/Cooling Rate 2 °C/min

Technical Parameters

Power Supply 90 - 230 V, 50 - 60 Hz, 100 W

LC Display target/actual temperature (simultaneously)

Temperature Adjustment with the included keyboard or via RS-232

Operation manually or software-controlled

Programming temperature gradient (99 steps or linear)

Leak Detector adjustable sensitivity; acoustic signal; turn-off switch

Safety self-check and auto-calibration at power-on; selectable turn-off temperature

Dimensions / Weight

150 x 470 x 300 mm (W x H x D) / 8.4 kg

Ordering Information

Order No.

A0585	Smartline Column Thermostat
A0527	Smartline Column Thermostat with 6-Port/2-Channel Injection Valve, 1/16", stainless steel
A1999	Smartline Column Thermostat with 6-Port/2-Channel Injection Valve, 1/16", PEEK

Technical data are subject to change without notice.

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Dr. Ing. Herbert Knauer GmbH