# WellChrom

## Filter-Photometer K-2001

## Filterphotometer K-2001

## Handbuch / User Manual

V7561, 06/2001

# CONTENTS

Using this Manual .....	4
Conventions in this manual.....	4
SOP's in this manual .....	5
The Filter-Photometer K-2001 .....	6
General Description .....	6
Optical Path of the Filter-Photometer K-2001 .....	7
Preparing the Filter-Photometer K-2001 for Operation .....	7
Unpacking.....	7
Standard delivery .....	7
Front Panel Controls .....	8
Front View of the Filter-Photometer K-2001 .....	8
Function of Foil Keys .....	8
Rear view .....	9
Power supply, ON/OFF, Selftest .....	9
Operating the Filter Photometer.....	10
Internal Software Structure .....	10
Installation of the Flow Cell.....	14
Capillary Connection to a HPLC System .....	14
Direct Control of the Filter-Photometer K-2001.....	15
Time programmed Chromatograms .....	16
Entering a Program.....	16
Running a Program .....	17
Connecting other Instruments to the Filter-Photometer K-2001.....	17
Using the remote control socket.....	17
Connections of the remote control socket .....	17
Assembling plug strips.....	19
Software Control of the Filter-Photometer K-2001.....	19
RS 232 Serial Interface .....	20
Simple Maintenance.....	21
Control of the lamp's functionality.....	21
Changing the lamp.....	21
Cleaning the flow cell .....	23
Analytical flow cells.....	23
Preparative flow cells.....	24
Adjusting the path length of the preparative flow cells.....	24
Flow cells with fiber optical connectors .....	25
Trouble shooting.....	26
Error messages and their reasons .....	26
Spare parts and accessories .....	27
Flow cells for the Filter-Photometer K-2001.....	27
Analytical Flow Cells.....	27
Preparative Flow Cells.....	27
U-Z View™ Micro Flow Cells .....	27
Spare Parts.....	27
Technical Data.....	28
Declaration of conformity.....	29
Guarantee statement.....	30

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## Using this Manual

This manual refers to the *WellChrom* Filter-Photometer K-2001 Firmware Revision 1.06 or higher. It is valid for any combination with analytical flow cells order number A 4061, A 4062, A 4063, A 4065, preparative flow cells order number A 4066, A 4067, A 4068, A 4069 and all UZ View™ micro flow cells in standard and fiber optic version.

### Conventions in this manual

Arrows like this: → ↔, used in block diagrams, indicate that the user is asked to press the corresponding arrow keys. The operation of arrow keys is defined as follows:

Cursor right: ▶      up: ▲      down: ▼      left: ◀.

**Important Hints** are marked by the marginal hand symbol.

**Special Warnings** are indicated by the marginal warning sign and printed in bold letters.

The marginal lamp symbol indicates helpful advice's.

## SOP's in this manual

The **Standard Operating Procedures (SOP)** provided with this manual offer a convenient way of structuring complex tasks in the operation of your Filter-Photometer K-2001. They include step-by-step instructions leading the user through all routine tasks during operation. They can be used for documentation purposes and be copied, applied signed, and filed in order to document the performance of the instrument.

**Please operate the instrument and all accessories according to instructions and SOP's in this manual. This ensures proper results and longevity of your equipment.**

<b>SOP 1</b>	Installation of the Flow Cell .....	14
<b>SOP 2</b>	Capillary connections .....	14
<b>SOP 3</b>	WAGO plug strip assembling.....	19
<b>SOP 4</b>	Changing the deuterium lamp. ....	22
<b>SOP 5</b>	Changing the halogen tungsten lamp.....	22
<b>SOP 6</b>	Purging the Flow Cell .....	23
<b>SOP 7</b>	Cleaning an analytical flow cell.....	23
<b>SOP 8</b>	Cleaning a preparative flow cell.....	24
<b>SOP 9</b>	Changing the path length.....	24

## The Filter-Photometer K-2001

### General Description

The *WellChrom* Filter-Photometer K-2001 is a valuable new detector, especially developed for routine measurements. The Filter-Photometer K-2001 satisfies all GLP requirements of a regulated laboratory.

The instrument is delivered in its standard version with a deuterium lamp as well as with 4 filters for the wavelengths 200, 220, 254, and 280 nm. Optional a tungsten-halogen lamp is available. Additional two further wavelength filters of your choice can be built in.

A rich palette of flow cells available for the KNAUER photometers, ranging from cells for Nano-HPLC with flow rates < 10 nL/min to preparative cells for up to 10 l/min, make the photometer K-2001 highly flexible in the full range of LC applications.

Moreover, a version of the Filter-Photometer K-2001 equipped with fiber optical connectors of the cell is available. This enables the spatial separation of detector and flow cell and thus the use of these instruments in hazardous locations.

After powering up the *WellChrom* Filter-Photometer K-2001 performs a self test automatically.

The high sensitive detector is marked by its excellent low noise ( $\pm 1 \times 10^{-5}$  AU) and also by a low baseline drift of  $5 \times 10^{-5}$  AU/h. The Autozero range is full scale.

The clear foil pad enables you to activate all basic functions in an easy to learn way. Furthermore you can enter, save, and run a time program for changing the filter. 10 program lines are available. The program can be executed as often as you want.

The control of the Filter-Photometer and the data collection occur digitally, to ensure a trouble free working of the instrument. Also an adjustable analog input and the option for an analog remote control are available.

### Optical Path of the Filter-Photometer K-2001

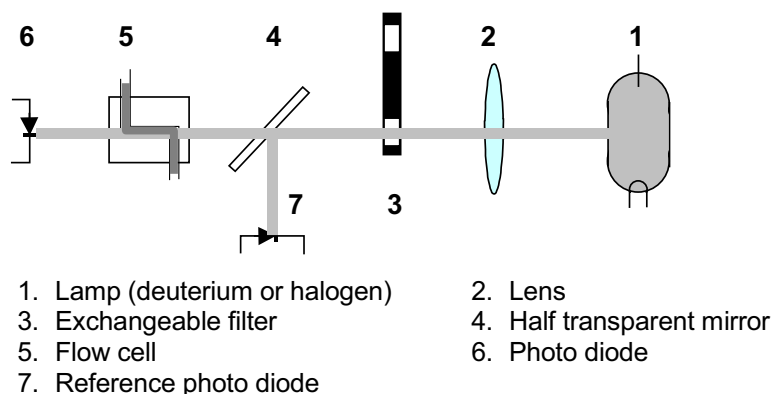
Fig. 1

*Optical Path of the Filter-Photometer K-2001*

The light emitted from the lamp (1) is focussed by the lens(2). The exchangeable filter (3) fades out the desired wavelength. The beam is then splitted by a half transparent mirror (4). One part beam delivers the reference signal (7) and the other after pathing the flow cell (5) the measurement signal (6).

## Preparing the Filter-Photometer K-2001 for Operation

### Unpacking

After unpacking, please check the device and accessories thoroughly for any damage that may have occurred during transport. If necessary, put forward any claim for damages to the carrier.

Use list "Standard delivery" and check that the Filter-Photometer K-2001 delivery is complete. Please contact our service department if you miss something or if you need support. Please fill out the guarantee registration card and return it to us immediately.

Remove the protection foils from the display and from the capillary inlet and outlet respectively.

### Standard delivery

WellChrom Spectro Photometer K-2501 without cell  
 Operation Manual  
 Power Supply Cable 230 V  
 RS-232 Cable  
 Integrator Cable  
 Plug Strips with Connectors  
 Analog Connection Cable

## Front Panel Controls

### Front View of the Filter-Photometer K-2001

1. Display information
2. Foil key area
3. Knurled cell screws
4. Flow cell housing
5. Flow cell cradle
6. Outlet
7. Flow cell
8. Path length of cell
9. Inlet

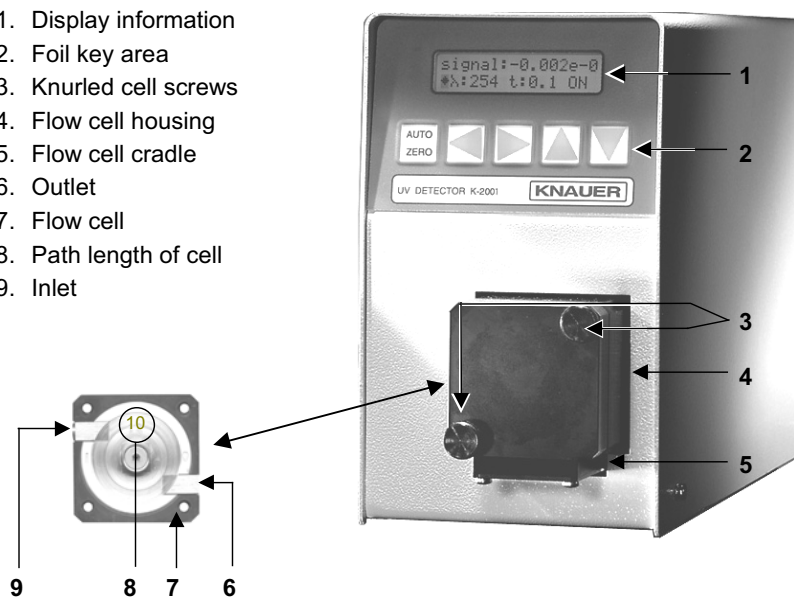
Fig. 2

Front Panel Elements of Photometer and Flow Cell

### Function of Foil Keys

The foil key area (2) in Fig. 2 „Front panel of the Filter-Photometer K-2001“ consists of an AUTOZERO key and four arrow keys.

#### AUTOZERO

Pushing this button will adjust the baseline. Usually the button should be pressed shortly before starting a chromatographic run. From any displayed menu the display returns to the SIGNAL-menu.

#### Arrow Keys

Use the yellow arrow keys  $\blacktriangleright$  „right“ or  $\blacktriangleleft$  „left“ to move and to position the cursor on the display as well as to confirm entered or selected values.

Using the yellow arrow keys  $\blacktriangle$  „up“ or  $\blacktriangledown$  „down“ you can change the selected parameter or the available options of it.

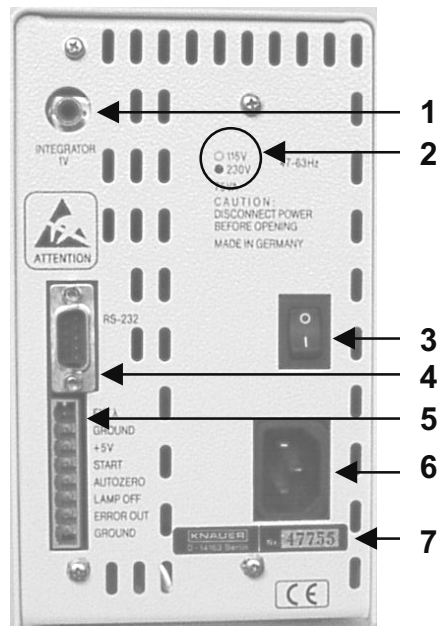
**Rear view**

Fig. 3

Rear view of the Filter-Photometer K-2001

- 1 Analog output (to recorder or integrator)
- 2 Indication of the input voltage (see Abb. 4)
- 3 ON/OFF switch
- 4 RS-232 Interface
- 5 Terminal strip Remote connections
- 6 Power connector
- 7 Serial number

**Power supply, ON/OFF, Selftest**

The Filter-Photometer K-2001 operates with 115 or 230 V AC of 47-63 Hz. The setting is done by the manufacturer on customers request. The standard setting is 230 V. If there is on any reasons the need of a change, please contact our service department. The actual setting is indicated on the rear panel of the instrument (Abb. 4).

Fig. 4 Indication of the input voltage

**Make certain that the correct voltage has been set on the rear panel of the instrument, the power supply is grounded and a corresponding 3-pole power cable is used.**

Connect the photometer to the power supply and switch on the instrument. The power switch is on the rear panel, see Abb. 3 on page 36.

After switching ON the display (item 1 in Fig. 2 on page 8) presents for a short time information's concerning the instrument version and number:

<b>*KNAUER</b>	<b>V 2.3</b>
<b>MINI-FIPHO</b>	

Any powering up the instrument includes a check of the electronics and an Autozero. Prior to this the lamp is heated to the working temperature for about 20 seconds indicated by the following display:

<b>signal:</b>	<b>t:0.2</b>	<b>-.----</b>
<b>◆λ:254</b>		<b>HEA</b>

The start routine will be finished by an automatically Autozero. Now the Filter-Photometer is ready to work, showing the display:

<b>signal:</b>	<b>t:0.2</b>	<b>0.0000</b>
<b>◆λ:254</b>		<b>ON</b>

## Operating the Filter Photometer

### Internal Software Structure

The software is divided in various menus, each of which allows particular settings and operational modes. You reach the single menu by positioning the cursor on the rhombus field **◆** followed by pressing the **Δ** or **∇** keys. As indicated in Abb. 5, the menus will be called up in an endless loop.

The single menu will be described in detail. Inside of any menu the cursor can be moved to the next or predecesing field using the **▷** or **◁** keys respectively. There it is possible to increase or decrease the corresponding parameter settings by help of the **Δ** or **∇** keys. In some cases you can scroll with the **Δ** or **∇** keys through the available options. Moving to an other entry field using the **▷** or **◁** keys respectively the entered changes will be confirmed.


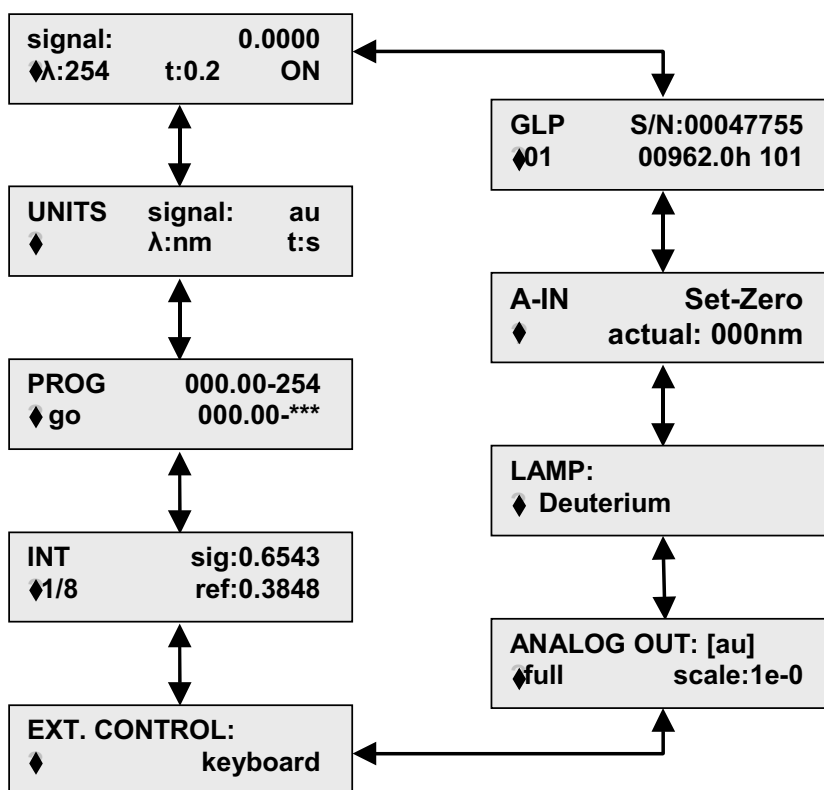
Fig. 5 Menu sequence of the Filter-Photometer K-2001

**The SIGNAL menu**

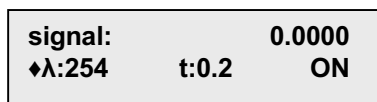The signal or main menu contains the data signal intensity **signal:**, wavelength **λ:**, time constant **t:** and the lamp status.

The actual measured signal value is shown in a fixed point manner with four digits. The measurement wavelength **λ:** can be selected out of the standard values 200, 220, 254, 280 nm, and if chosen of your optional filters.

Using a time constant you can achieve a signal smoothing. The time constant **t:** can be set to the values of **0.1, 0.2, 0.5, 1, 2, 5** or **10** seconds. The larger this value is set the more the signal will be smoothed. A time constant of 1 s fits most of analytical purposes best.

At last you can switch ON/OFF the lamp in this menu, indicated by **ON** and **OFF** respectively.

**The UNITS menu**

This menu is for information only, no changes are possible. The units of the signal menu values are shown.

### The PROG menu

<b>PROG</b>	<b>000.00-254</b>
◆ <b>go</b>	<b>000.00-***</b>

The PROG menu serves to enter a time based program for changing the wavelength's. The way how a program is entered and runned will be described in the section Entering a Program on page 16. Inclusive both starting and ending lines a program can consists of 11 lines maximum.

### The INT menu

This menu is for information only, no changes are possible.

<b>INT</b>	<b>sig:0.6543</b>
◆ <b>1/8</b>	<b>ref:0.3848</b>

The values **sig** and **ref** record the light intensities of the signal and reference channel respectively. The **sig** and **ref** values are in the range between 0 and 1. They depend on the selected wavelength, the solvent in use, the flow cell as well as on the age of the lamp. The ratio of both the values should be between 1:1.5 and 1:1.8. Significant deviations are an indicator for instance for a dirty lens or a to strong absorption of the solvent at the selected wavelength.

The value **1/8** behind the rhombus in the second line is a measure for the integration time and it is only of interest in case of servicing.

### The EXT. CONTROL menu

<b>EXT. CONTROL:</b>	
◆	<b>keyboard</b>

This menu enables you to select the control mode: either from other instruments (external control = **analog**) or in a **stand-alone** mode (selection **keyboard**)

### The ANALOG OUT menu

<b>ANALOG OUT:</b>	<b>[au]</b>
◆ <b>full scale:</b>	<b>1e-0</b>

The ANALOG OUT menu is used for calibration of the analog output signals. Generally the maximum output signal is 1 V. In accordance with your actual conditions you can determine how many AU correspond to this signal maximum or **full scale** deflection. The values of 1e-4, 2e-4, 5e-4, 1e-3, 2e-3, 5e-3, 1e-2, 2e-2, 5e-2, 1e-1, 2e-1, 5e-1, 1e-0, 2e-0, 5e-0 or 1e+1 AU (corresponding to 0.0001, 0.0002, 0.0005, 0.001, 0.002, 0.005, 0.01, 0.02, 0.05, 0.1, 0.2, 0.5 1, 2, 5 and 10 AU respectively) are selectable.

### The Lamp menu

This menu is for entering the type of lamp used to adapt the software. The options Deuterium and Halogen are possible.

LAMP: ◆ Deuterium
----------------------

The setting shall only be changed, if really a change of the lamp type was carried out. If the setting and the actual installed lamp type are not in agreement, the instrument will not work properly.

### The A-IN menu

A-IN	Set-Zero
◆ actual:	000nm

The A-IN-(Analog-IN) menu is used to calibrate the analog input for automatic adjusting the monochromator. Activating the **Set-Zero** entry function interprets a voltage signal at the input as 000 nm. If then a signal of for instance 5 V is applied to the input, the **actual** field will show 500 nm as recognized wavelength value, considering a setting of 1 V = 100 nm, 500 nm. To produce a second calibration point the displayed wavelength can be changed pressing the ▽ or Δ keys.

### The GLP menu

GLP	S/N:00047755
◆01	00962.0h 101

The GLP menu contains the following information: serial number total number of lamps used, total operating hours of the current lamp and number of lamp starts.

Pressing the ▷ or ◀ keys you access the lamp counter. Pressing now one of the Δ or ▽ keys, the lamp counter will be **incremented** in both cases by 1. At the same time the ignition counter as well as the operating time counter will be set to zero!

This entry is irreversible! Make certain, that this function is used only after the lamp was really exchanged.

## Installation of the Flow Cell

The Spectro Photometer K-2501 comes factory configured without a flow cell. The device is equipped with a "dummy" cell which does not have any optical parts. Before operating the photometer it is necessary to install an appropriate KNAUER flow cell.

The instrument or the lamp are not needed to be switched off for changing the flow cell. When the flow cell is removed an error message **intens.overflow** will appear in the upper display line because the analog to digital converter gets an overflow due to the incoming light. This is **not** an instrumentation **fault!** After having installed the new flow cell properly, press any button to delete the error message.

### SOP 1

#### Installation of the Flow Cell

This instruction refers to the Filter-Photometer K-2001 without fiber optical connectors. Compare Fig. 2 „Front Panel Elements of Photometer and Flow Cell“ on page 8.

Loosen the two knurled cell screws (3) by hand and remove them.

1. Pull out the flow cell housing (4).
2. Take the cell or dummy cell with two fingers and remove it upward.
3. Insert the new flow cell (7) and make sure as well that the engraved specification (8) points towards the user and can be read as the fixing hole on the back side of the cell meets the corresponding metal pin of the photometer's housing.
4. Now push the complete system towards the housing, insert the two screws and tighten them manually.

## Capillary Connection to a HPLC System

**Before taking a measurement cell filled with fluid into operation, please make certain that the used eluent is miscible with that one used previously. Otherwise purge the flow cell with a medium miscible with both the other fluids.**

### SOP 2

#### Capillary connections.

This instruction refers to the Filter-Photometer K-2001. Compare Fig. 2 „Front Panel Elements of Photometer and Flow Cell“ on page 8.

5. Connect the outlet of the HPLC column to the inlet bushing of the flow cell (9).

Please use DYNASEAL bushings and the shortest possible capillary with small internal diameter in order to keep the dead volume as small as possible.

6. Push the bushing, the clamping ring, and the sealing ring onto the capillary. Please take care on the sequence and orientation of fittings, see Fig. 6 „DYNASEAL Capillary connections“.
7. Push the capillary as far as possible into the flow cell input.
8. Fasten the bushing by hand.
9. Connect the flow cell outlet (6) using a capillary or teflon tube (ID > 0,5 mm) to a waste bottle.

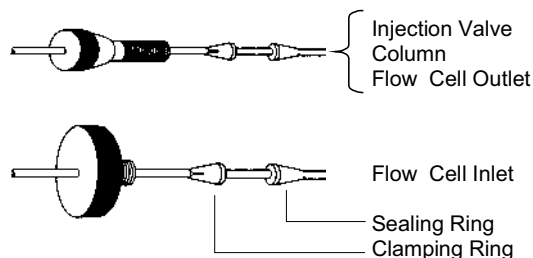
Fig. 6 DYNASEAL Capillary connections

The capillary in a simple isocratic system is shown in the following figure. For perspective reasons the Filter-Photometer K-2001 is symbolized by the flow cell and the used HPLC pump by its pump head.

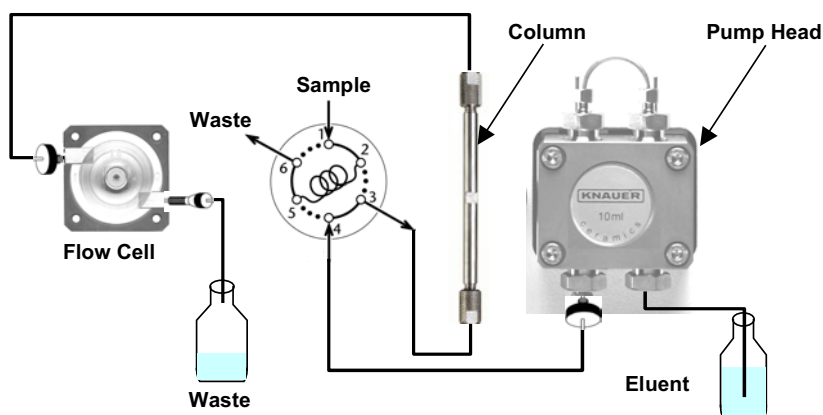
Fig. 7 Isocratic HPLC System

### Direct Control of the Filter-Photometer K-2001

In this section it will be described, how the Filter-Photometer K-2001 is to configure and to operate.

Switch the instrument on keeping in mind the hints given in the section Power supply, ON/OFF, Selftest on page 9.

**Prior to the first measurement wait for about 15 minutes for warming up the instrument with the HPLC pump also switched on. In case of especially sensitive measurements even prolonging this warming up time may be needed.**

If the instrument after passing the start routine is displaying

**Instrument is in EXTERNAL mode!**

it is under extern control by other devices. To control the photometer directly please change the setting in the EXT. CONTROL menu (see page 12) to **keyboard** control.

The lamp starts generally with any powering up automatically. However at any time it can be switched of and on in the SIGNAL menu (page 11).

Also in the SIGNAL menu select the desired wavelength and the appropriated time constant. In principle your Filter-Photometer K-2001 is now ready to take simple chromatograms.

## Time programmed Chromatograms

### Entering a Program

Programs are used to set and to change the wavelength during a run. Activate as described on page 12 the PROG menu. If no program is already saved you will get the following display:

<b>PROG</b>	<b>000.00-254</b>
<b>◆ go</b>	<b>000.00-***</b>

In the first line of the program you always will find **000.00** (0 minutes, 0 seconds) as starting point and an actual wavelength. The proceeding line contains the same time and "\*\*\*\*" for the wavelength. This ending line is used to create new lines for the program.

The cursor can be moved to the next or predecessing field using the  $\triangleright$  or  $\triangleleft$  keys respectively. There it is possible to increase or decrease the corresponding parameter settings by help of the  $\Delta$  or  $\nabla$  keys. Moving to an other entry field using the  $\triangleright$  or  $\triangleleft$  keys respectively the entered changes will be confirmed.

**You can leave the entry fields without confirming the changes by pressing the AUTO-ZERO button. The cursor moves back to the menu selector (rhombus field).**

First enter a new time in the editing line. The wavelength of the predecessing line will be adopted for possible changes. Only than the wavelength field of this line is accessible for changes. Simultaneously a new ending line is created, which can be achieved by moving the cursor with the  $\triangleright$  key. It is impossible to enter a time value smaller than the previous one. This ensures programs without time conflicts.

Including both the start and the ending line a program can consists of 11 lines as a maximum. Therefore you can measure during a single run with up to 10 different wavelength's.

**Caution! If you enter in any existing program line the time value of the predecessing line it becomes automatically the ending line. That means all following lines will be irreversible deleted!**

To delete a program enter in its second line the time value **000.00** (0 minutes, 0 seconds).

## Running a Program

A program can be executed via an external signal (section *The PROG menu* on page 12), as well as manually. Position the cursor on the **go** field and press one of the  $\Delta$  or  $\nabla$  keys.

A running program is indicated by **RUN**, see the following display, accompanied by the actually measured **signal**, the running time and the actual wavelength.

signal:	0.4567
RUN	001.30-254

Any programmed change of the wavelength is followed by an AUTO-ZERO.

A running program can be stopped at any time by pressing of any key. If you leave a program the last wavelength setting will remain active. A program will not be lost by switching off the instrument.

## Connecting other Instruments to the Filter-Photometer K-2001

### Using the remote control socket

On the rear panel side of the Filter-Photometer K-2001 an electrical connector socket is located (item 5 in Fig. 3 on page 9) which serve to send or receive signals from other instruments. For example outgoing start signals of an injection valve or an autosampler can be laid on the START input. All voltages have to be mounted between GROUND and the corresponding event.

**Please avoid touching the electrical contacts of the socket lines. Electrostatic discharges when touching the contacts could damage the electronics of the device.**

To operate the photometer external controlled change the setting in the EXT. CONTROL menu to **analog** (see page 12).

### Connections of the remote control socket

Two of the eight positions on the remote control socket are **ground** connections one is for ERROR OUT and four serve as control connections.

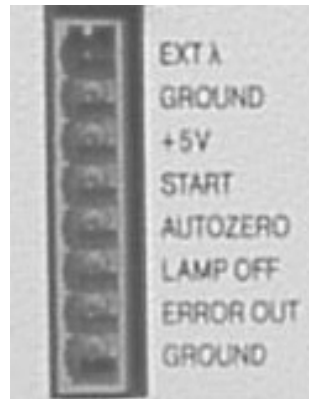
Fig. 8

## Remote Control Connections

- EXT λ** Control voltage signal for automatic wavelength selection (maximum voltage: 10 V = 10 mV/nm).
- +5V** **Caution! This connection is not allowed to use. It serves for service purposes only.**
- START** Short circuit to GROUND causes a program execution (see *Running a Program* on page 17).
- AUTOZERO** Short circuit to GROUND triggers an Autozero signal. Measurement restarts after the signal is switched off.
- LAMP OFF** Short circuit to GROUND means lamp **OFF**. Open input means lamp **ON**.
- ERROR OUT** Error signal (TTL, open collector), remains active as long as an error will be displayed e.g. the lamp does not start.

The Fig. 9 shows as an example the wiring of the Filter-Photometer K-2001 with the KNAUER Interface box.

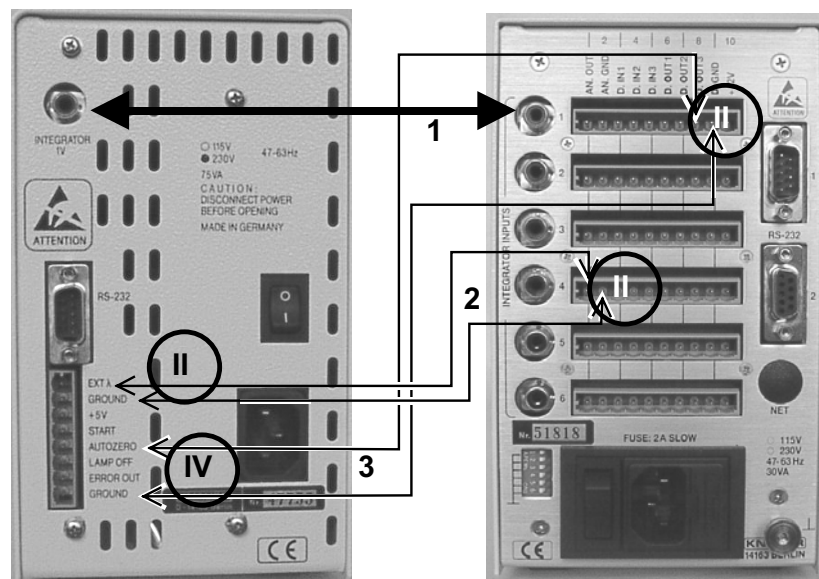
Fig. 9

## Connections of the Filter-Photometer K-2001 with a KNAUER-Interface box

The connection **1** (Integrator output) is mounted using a cinch-cinch-connection cable. The connections **2** and **3** are closed with cables you are obliged to mount yourself as described in the following section. The connection **2** needs two 2-pole plugs (II) and the connection **3** needs one 2-pole plug (II) **and** one 4-pole plug (IV).

### Assembling plug strips

For wiring the Filter-Photometer K-2001 with other instruments you need cables with WAGO plug strips. Plug strips with 2, 3 or 4 positions are included in the standard delivery. They are mounted as follows:

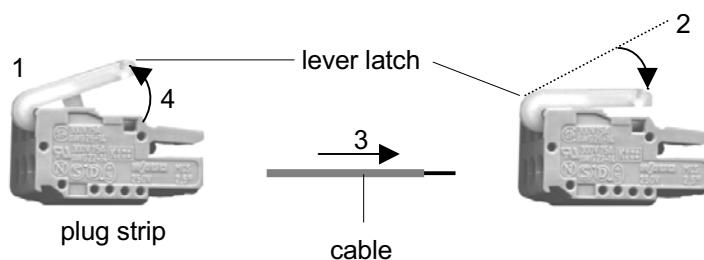
Fig. 10

Assembling plug strips

### SOP 3

#### WAGO plug strip assembling

1. Insert the rounded end of the lever latch into the square opening of the selected connector of the plug strip.
2. Press the catch down as indicated by arrow.
3. Insert the uninsulated end of the cable into the opening under the catch.
4. Release the catch and remove the lever latch from the plug.

The cable is now firmly anchored in the plug strip.

### Software Control of the Filter-Photometer K-2001

The full capabilities of the Filter-Photometer K-2001 are accessed under operation with the HPLC software packages EuroChrom<sup>®</sup> for Windows or ChromGate<sup>®</sup>.

Fig. 11

HPLC software packages

This chapter gives only a short information regarding the connections of the detector when working with the Filter-Photometer K-2001 under EuroChrom<sup>®</sup> or ChromGate<sup>®</sup> software control. For more detailed information's concerning the features of the software, please consider the users handbook of the software.

**Working under software control, the PROG-menu must be empty, no program is allowed!** Only then a START signal will be interpreted as command to start the measurement. In case a program is entered, any START signal will be interpreted as command to start this program.

### RS 232 Serial Interface

The RS 232 serial interface on the rear side of the device, item **4** in Fig. 3 on page 9, enable digital data transfer between the Filter-Photometer K-2001 and a PC, equipped with HPLC software (EuroChrom<sup>®</sup> or ChromGate<sup>®</sup>). Please connect this interface directly or if necessary via an interface multiplier to the **COM-Port** of your computer.

To operate the photometer external controlled by a software package change the setting in the EXT. CONTROL menu to **keyboard** (see page 12).

## Simple Maintenance

### Control of the lamp's functionality

The deuterium lamp used with the photometer K-2001 has an extended life time to ensure long-time functionality and reliable measurements with low noise and baseline drift as well as high sensitivity. The actual using time of the lamp depends on different factors, like the number of lamp starts, the average burning time and your requirements concerning noise and sensitivity.

To check the functionality of the lamp, the two intensity values **sig** and **ref** to be found in the SIGNAL menu on page 11 provide helpful information. The ref value refers to the intensity of the light measured in the reference channel and can be used for checking the quality of the lamp.

On delivery of the photometer, this value is set to the range of 0.39 to 0.78, measured with the dummy cell at a wavelength of 254 nm.

We recommend to check the **ref** value at regular intervals under the conditions mentioned above (dummy cell,  $\lambda = 254 \text{ nm}$ ). This applies especially in case higher noise levels or decreased sensitivity are observed on working with the photometer K-2001. In these observations coincide with a **ref** value of approx. 0.1 or less, a new deuterium lamp should be installed.

### Changing the lamp

**Remove the power plug before opening. Please let the lamp cool down for at least 15 minutes after switching it off.**

To change the lamp see Fig. 12 to Fig. 14.

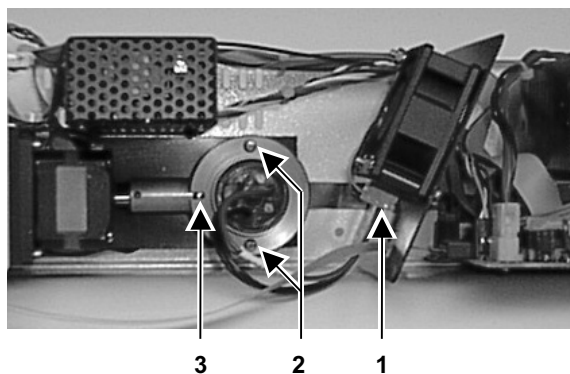
Fig. 12

Top view of the Filter-Photometer K-2001 with removed housing

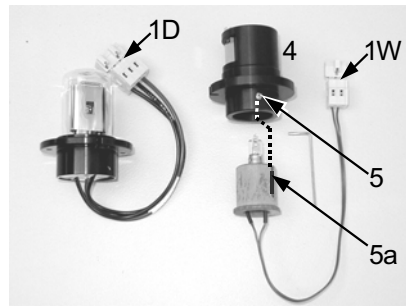
Fig. 13 Lamp assembly

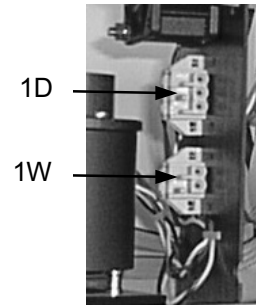
Fig. 14 Lamp sockets

**Do not touch the glass of the lamp. Should you touch it accidentally, clean it thoroughly with a lint free cloth and i-propanole.**

#### SOP 4

Changing the deuterium lamp.

This instruction applies to the Filter-Photometer K-2001.

1. Unscrew the housing and remove it by lifting.
2. Pull the 3-pole plug of the old lamp (item 1D). Unscrew the two screws in the lamp socket (item 2) with a screw driver and remove the whole lamp from the instrument including its cables.
3. Insert the new lamp assuring that it is correctly seated in the guiding slot (item 3).
4. Screw it in securely (item 2) and connect the 3-pole plug into the 3-pole socket (item 1D).
5. Increase the lamp counter in the GLP-menu. The operating time counter will be set to zero by this implementation.
6. Put the lid back on from the top of the instrument and fasten the screws of the cover.
7. Check the intensity values in the INT-menu.

**Hint: If a new lamp is installed, it takes about 24 hours to reach the optimum working conditions.**

#### SOP 5

Changing the halogen tungsten lamp.

This instruction applies to the Filter-Photometer K-2001.

1. Unscrew the housing and remove it by lifting.
2. Pull the 2-pole plug of the old lamp (item 1T). Unscrew the two screws in the lamp socket (item 2) with a screw driver and remove the whole lamp from the instrument including its socket (item 4) and cables.
3. Loosen the pin screw (item 5) and pull the lamp out of its socket.
4. Insert the new lamp assuring that it is correctly seated in the guiding slot (item 5, 5a) and fasten the pin screw.
5. Insert the socket with lamp into the instrument, screw it in securely (item 2), and connect the 2-pole plug into the 2-pole socket (item 1T).
6. Increase the lamp counter in the GLP-menu. The operating time counter will be set to zero by this implementation.

7. Put the lid back on from the top of the instrument and fasten the screws of the cover.
8. Check the intensity values in the INT-menu.

### Cleaning the flow cell

Noisy baselines and low sensitivities may be due to a dirty flow cell. This may also be indicated by a low value for **sig** in the signal menu when flushing the cell with pure solvent. In most cases it is sufficient to purge the flow cell according to the following SOP.

#### SOP 6

##### Purging the Flow Cell

1. Purge the flow cell using one of following solvents: sodium dodecyle sulfate (SDS), 1m HCl, 1m NaOH, ethanol, or acetone.
2. Run the solvent through the flow cell using a syringe and leave for approximately 5 minutes..
3. Rinse extensively with water and the blow dry using a gentle stream of pure nitrogen.

**Never dry with compressed air from a „house“ line as this will contain microdroplets of oil that will coat the cell.**

When the optics module is not in use, disconnect the flow cell and clean out traces of salt and protein with a syringe filled with distilled water. Before storing the flow cell inject a dilute solution (10-25%) of ethanol or i-propanol to prevent microbial growth.

In case the flow cell purging do not provide sufficient success, all flow cells can easily be disassembled for cleaning the lenses.

### Analytical flow cells

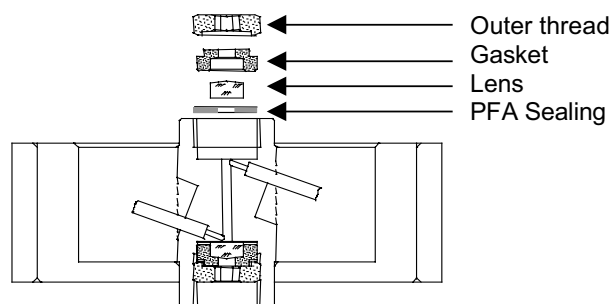
Fig. 15

Sectional view of an analytical flow cell

#### SOP 7

Cleaning an analytical flow cell. This instruction applies to the analytical flow cells A4061, A4062, A4063, and A4065.

1. Unscrew the outer threads with the 3 mm hexagonal spanner enclosed in the flow cell's delivery.
2. Remove the black gasket that carries the lenses with a pair of tweezers or by gently tapping it on a clean surface. The lens is embedded in the gasket and sealed against the flow path with a PTFE seal. This seal should be changed every time when disassembling the flow cell.
3. Take out the lenses and clean them by wiping them with a soft cloth or with an appropriate solvent in an ultrasonic bath. Be careful not to touch the clean lenses with the fingers.

4. Reassemble the cell in the reverse manner, making sure that the PTFE seal does not block the light path.
5. Tighten the outer threads carefully with the spanner in order not to damage the lenses.

### Preparative flow cells

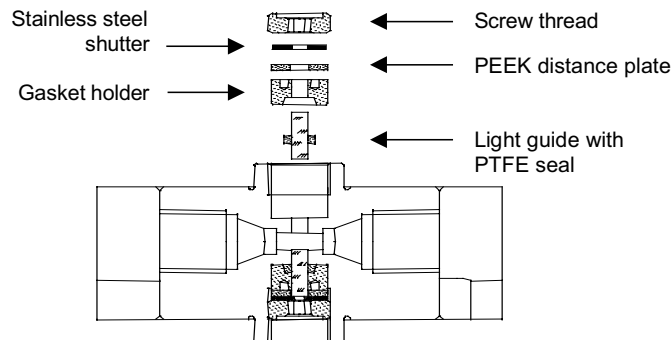
Fig. 16 Sectional view of a preparative flow cell

**SOP 8** Cleaning a preparative flow cell. This instruction applies to the preparative flow cells A4066, A4067, A4068, and A4069.

The preparative flow cells have a rod shaped light guide instead of the concave lens in the analytical cells.

1. Unscrew the outer thread with a hexagonal spanner
2. Take out the stainless steel shutter and the PEEK distance plate (not present in A4069).
3. Get hold of the gasket holder with the light guide by grasping it with a pair of tweezers, using the indentations on the outer side of the gasket holder.
4. Push out the light guide and strip the PTFE sealing ring in order to clean the lens.
5. Reassemble the cell in reverse order. Use a new PTFE sealing ring after every disassembly to ensure the consistence of the flow cell.

### Adjusting the path length of the preparative flow cells

**SOP 9**

Changing the path length.

This instruction applies to flow cells A4066, A4067 and A4068 with 1/8" and 1/4" connectors. Path lengths can be adjusted to 2, 1.25 and 0.5 mm. On delivery the path length is set to 2 mm. To reduce the path length to 1.25 or 0.5 mm, follow the instructions given.

1. Unscrew the outer thread with the 3 mm hexagonal spanner.
2. Take out the stainless steel shutter and the PEEK distance plate.
3. Remove the distance plate, put back the stainless steel shutter and refasten the thread carefully.

With the distance plate removed, the rod shaped light guide is pushed further into the flow cell (0.75 mm), thus resulting in a reduced path length of 1.25 mm. To reduce the path length further to 0.5 mm, follow the same procedure on the other side of the cell.

To extend the path length again in steps of 0.75 mm, the PEEK distance plates have to be inserted again.

4. Loosen the outer thread, remove the stainless steel shutter and take out the gasket holder by using a pair of tweezers.
5. Push the light guide approximately 1 mm to the outside to enlarge the path length. Use a clean cloth and do not touch the light guide with the fingers.
6. Put the gasket holder back into the cell.
7. Insert the PEEK distance plate again and then the shutter.
8. Fasten the outer thread carefully.

When fastening the threads, the rod shaped light guide is pushed back in the correct position inside the cell. Inserting a distance plate thus enlarges the path length for 0.75 mm. It is not necessary to change the PTFE sealing ring when adjusting the path length.

### Flow cells with fiber optical connectors

Changing the path length of flow cells with fiber optical connectors can be done in the same way. In these cells the screwthread is replaced by a special adapter (Fig. 17). They additionally contain a lens to focus the light onto the optical fibers.

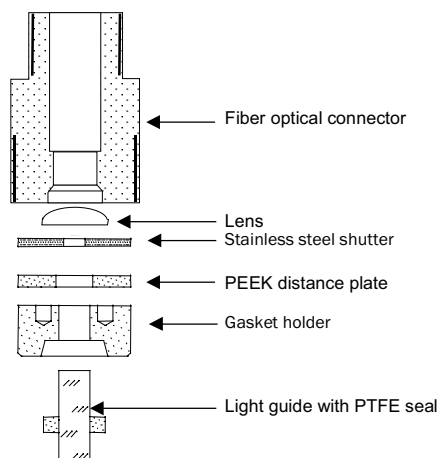
Fig. 17

*Fiber optical connector of a preparative flow cell*

## Trouble shooting

### Error messages and their reasons

Error message	probable cause	solution
<b>Deuterium lamp don't start</b>	Lamp defect	Exchange lamp
	Halogen lamp installed with lamp menu setting <b>Deuterium</b>	Change the setting in lamp menu
<b>ERROR zero position!</b>	Zero position can not be found	To perform a calibration switch the instrument OFF and ON
	Deuterium lamp installed with lamp menu setting <b>Halogen</b>	Change the setting in lamp menu
<b>Intensity overflow</b>	Light intensity too high, e.g. during exchanging the flow cell	Push any key after proper installing the flow cell
<b>ERROR: moving filter wheel</b>	The filter wheel is blocked for any reason	If this reason is not obviously, contact our service

If only **ERROR** is displayed press any key to continue. If there occur repeated problems starting the lamp or with calibrations please contact our service department.

The message **ERROR** interrupts a running program of your Filter-Photometer K-2001 as long as you have not pressed any key or an external start signal executes the program newly. The message **signal overflow** do not cause a program interruption.

## Spare parts and accessories

### Flow cells for the Filter-Photometer K-2001

All flow cells are also available equipped with fiber optical connectors for the use with the fiber optics version of the photometer K-2001.

#### Analytical Flow Cells

Order No. Cell type	Layer Thickness (mm); Connector	ID channel (mm)	Volume ( $\mu$ L)	Material	Flow Range (mL/min)	Maximum Pressure (bar)
A4061	10 mm; 1/16"	1,1	10	stainless steel, with heat exchanger	20	300
A4042	3 mm; 1/16"	1,0	2	stainless steel	50	300
A4045	3 mm; 1/16"	1,0	2	PEEK	50	30

#### Preparative Flow Cells

A4066	0,5/1,25/2 mm 1/8"			stainless steel	1.000	200
A4067	0,5/1,25/2 mm 1/8"			PEEK	1.000	100
A4068	0,5/1,25/2 mm 1/4"			stainless steel	10.000	200
A4069	0,5 mm 1/16"			stainless steel	250	200
A4095	0,5 mm 1/16"			PEEK	250	100

#### U-Z View™ Micro Flow Cells

A4091	8 mm 1/16"	0,150	0,140	fused silica	0,10	500
A4092	8 mm 280 $\mu$ m	0,015	0,035	fused silica	0,01	500
A4093	8 mm 280 $\mu$ m	0,020	0,003	fused silica	0,001	500
CE Cell: A4097	1 mm 280 $\mu$ m			stainless steel		

#### Spare Parts

A4071	Deuterium lamp
A4072	Halogen lamp
A4073	Halogen lamp with socket (needed for first installation of a halogen lamp)
M1642	Power supply cable
A0884	RS-232 connection cable
A1402	Set of plug strips (3x4; 2x3; 1x2 connections) including lever latch
G1023	Integrator cable
M1588	Analog connector cabal
A1467	10 pin Ribbon Cable
A1131	Repair kit for analytical flow cells
A1132	Repair kit for preparative flow cells

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### Manufacturer's address:

Hegauer Weg 38  
14163 Berlin, Deutschland

**WellChrom Filter-Photometer K-2001**, Order Number **A 4160** and  
**A 4200** (fiber optical version)

complies with the following requirements and product specifications:

- Low Voltage Ordinance (73/23/EWG);  
EN 61010 – 1 (1993)
- Engineering Guidelines (89/392/EWG)
- EMV Ordinance (89/336/EWG)  
EN 50081 – 1 (1992)  
EN 55011 (1991) Class B  
EN 55022 (1987) Class B  
EN 50082 – 1 (1992)  
IEC 801 – 2 (1984),  
includes IEC 41 B (sec) 81 (1992)  
IEC 801 – 3 (1984)  
IEC 801 – 4 (1988)

The product was tested in a typical configuration.

Berlin April, 10<sup>th</sup> 2000

Bernward Rittgerodt (Managing Director)

The CE Shield is attached to the rear of the instrument.

## Guarantee statement

The guarantee period of the WellChrom Filter-Photometer K-2001 is 12 months beginning from the date of dispatch from Berlin. Operation inconsistent with manufacturer's instructions or damage caused by unauthorized service personnel are excluded from guarantee. Damage caused by blockages and wear and tear parts such as fuses and seals are not covered by the guarantee. Claims under this guarantee are valid only if the enclosed guarantee card is returned to us at the address shown below within 14 days of receipt of the instrument. Defective pumps should be sent to the manufacturer for repair.

Wissenschaftliche Gerätebau  
Dr. Ing. Herbert Knauer GmbH  
Hegauer Weg 38  
D-14163 Berlin  
Tel: 030 – 809 727 – 0  
Fax: 030 – 801 50 10  
e-Mail: [info@knauer.net](mailto:info@knauer.net)  
[www.knauer.net](http://www.knauer.net)

If we find a defect covered by the guarantee, repair or replacement, at our discretion, will be carried out free of charge. Packing and transport costs are borne by the purchaser.

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