DAD K-2800

First Steps Erste Schritte

V7601, 01/2003

Version 1.0

Contents

| | |
|--|----|
| Introduction | 3 |
| Delivery Contents | 3 |
| General description | 4 |
| How to install DAD K-2800 | 6 |
| PC Specifications | 6 |
| Installation of the DAD K-2800 | 7 |
| Operating the Detector with ChromGate | 9 |
| Detector setup | 9 |
| Settings and initial tests | 10 |
| Operating the Detector with EuroChrom, Preparative Version | 13 |
| Detector setup | 13 |
| Settings and initial tests | 15 |
| Spare Parts and Accessories: | 17 |
| Technical Data | 17 |
| Declaration of conformity | 18 |
| Guarantee statement | 19 |

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Introduction

The KNAUER Diode Array Detector K-2800 is a top performance UV/VIS/NIR-detector based on highly advanced fiber optical technology. It features very wide spectral range, high sensitivity, very low noise level, and very low drift.

There are two versions of the detector: DAD K-2800 UV/VIS/NIR (A4444) the full wavelength range diode array detector with D₂-lamp and halogen lamp installed and the DAD K-2800 UV (A4445) with only a D₂-lamp installed (190-740 nm range). The DAD K-2800 UV can be easily upgraded to the DAD K-2800 UV/VIS/NIR by authorized KNAUER service engineers.

Due to its unique design the KNAUER Diode Array Detector K-2800 requires no moving parts, no reference channel, no special 'see through' lamps and no *power on*-calibration. It is a fully software controlled detector which means that all functions and settings of the detector are directly accessible via KNAUER ChromGate or EuroChrom Preparative Version.

The whole range of KNAUER flow cells from analytical cells (e.g. 1/16", 10 mm, 10 µl, 20 ml/min, 300 bar) up to large preparative cells (e.g. 1/4", 1/1.25/2 mm, 10,000 ml/min, 200 bar) is available for the DAD K-2800. By simply connecting longer fiber optical cables, the flow cell can be moved from inside the instrument to any other place – for example for explosion proof applications.

The KNAUER Diode Array Detector K-2800 can be used to record up to four chromatograms as well as up to 10 spectra per second (scan range can be chosen within 190-1024 nm). Together with the chromatography software DAD K-2800 provides 2D and 3D chromatograms, spectral library and peak purity functionality. The high amount of data which is sent to the PC is handled by a special high speed RS 422 interface card. For preparative applications the KNAUER Diode Array Detector K-2800 offers two adjustable analog outputs.

Delivery Contents

The KNAUER Diode Array Detector K-2800 is delivered with:

- PCI-card for PC/DAD communication
- RS-422 serial communication cable to connect PCI-card and DAD
- Mains cable

Flow cells can be chosen from the wide range of analytical and preparative KNAUER flow cells. See Spare Parts and Accessories on page 17.

General description

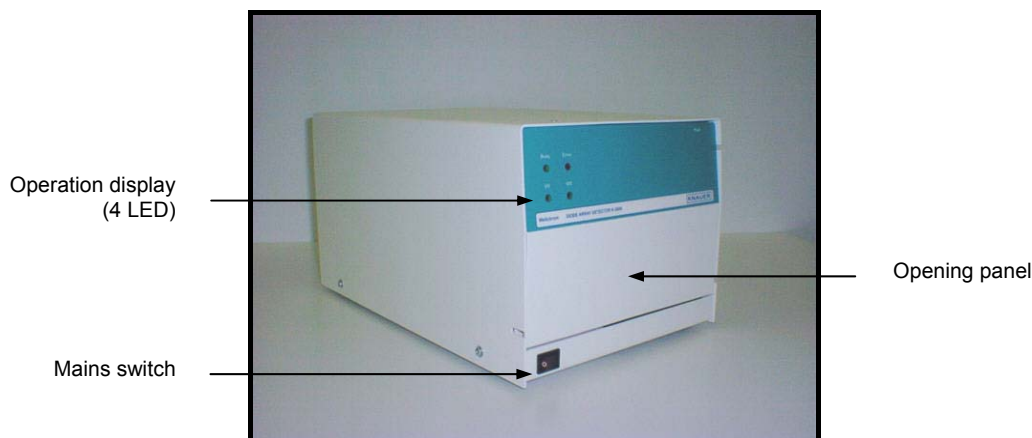
Fig. 1 Front panel view

The DAD K-2800 is a fully software controlled instrument which needs only the mains switch and four LEDs on the front.

When the DAD K-2800 is switched on the integrated light in the mains switch will turn on and the red LED in the operation display will flash once after about two seconds. The green **UV-**, **VIS-** and **Busy-**LEDs will only light up when the detector is connected with a PC and initialized by the chromatography software.

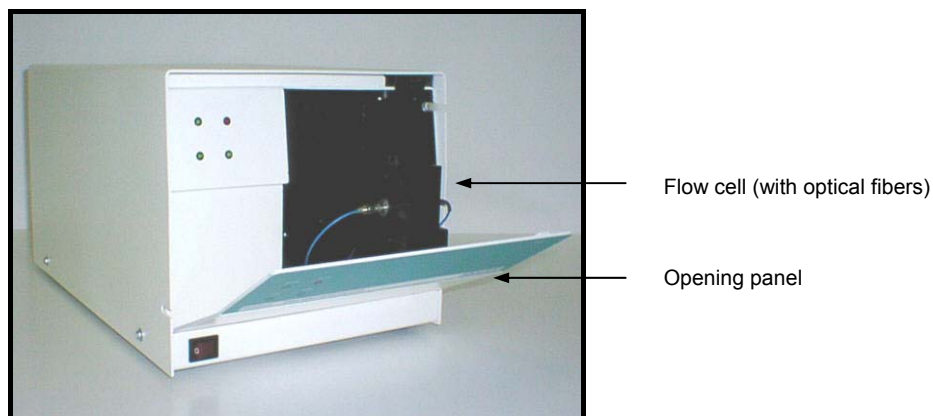
Fig. 2 Front panel view (open, with cell)

The flow cell, capillaries and fiber optical cables are located behind the front panel. The flow cell is carried by the moveable flow cell holder. For best results, the front panel should be closed during operation of the detector.

The flow cell is put into the flow cell holder which can be pulled out for easy access to the flow cell. Flow cells should be installed with capillary inlet down and capillary outlet up to prevent bubbles in the flow cell. (KNAUER analytical flow cells are marked with 'I' at the inlet and 'O' at the outlet. For this type of cells inlet and outlet can not be exchanged because this would result in increased noise. For all other cells inlet and outlet are identical.) It is recommended the inlet capillary be installed after the cell is placed in the cell holder. For preparative (1/8") cells specially prepared connecting capillaries are supplied with the cells. 1/4" cells can not be placed in the cell holder.

The detector's two fiber optical cables are connected with the cell's connectors and are screwed tightly without using any tool. The fiber optical cables should be arranged in a way that they are bent as least as possible.

Fiber optical cables are very sensitive to bending and may break easily. They must be handled very carefully and may only be bent as much as necessary for connecting the flow cell.

A broken fiber optical cable causes very low light intensity which leads to very high noise and drift.

As the life time of fiber optical cables depends very much on it's handling by the user, these cables, as well as lamps, are normally not covered by any warranty (except the user can prove that it is not his fault or that the instrument was already delivered with a bad fiber optical cable or lamp).

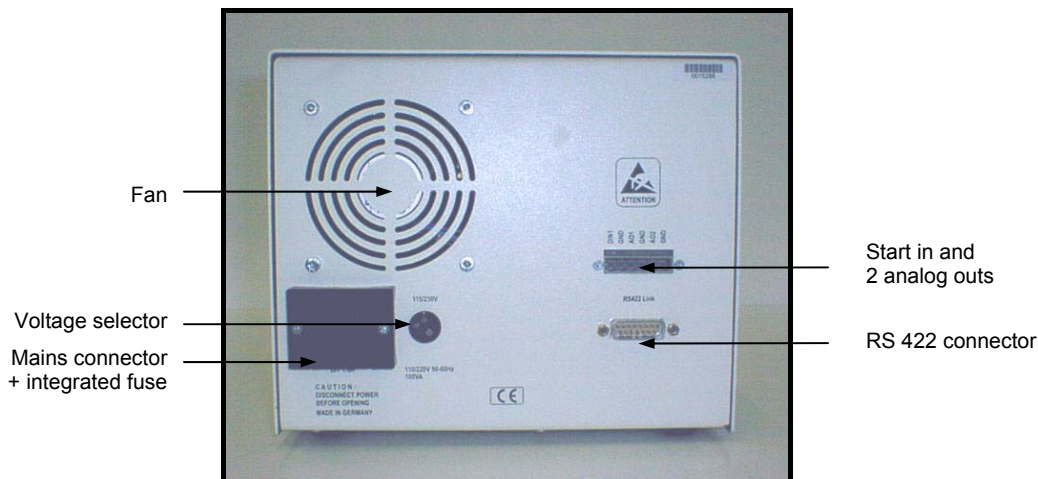
Fig. 3 Rear panel view

All available electrical connectors are located on the instrument's rear side. The RS-422 socket is needed to connect the DAD K-2800 to a PC. Start in is used to send a start signal to the detector (trigger must be set to external and DAD-K-2800). The two scalable analog outputs are only activated in combination with EuroChrom Preparative Version. Analog out 1 represents channel 1 and analog out 2 represents channel 2. The voltage selector sets the instrument either to 220-230 V or 110-115 V. Default setting is 220-230 V.

Before using the instrument please check whether the setting of the voltage selector fits with your local voltage.

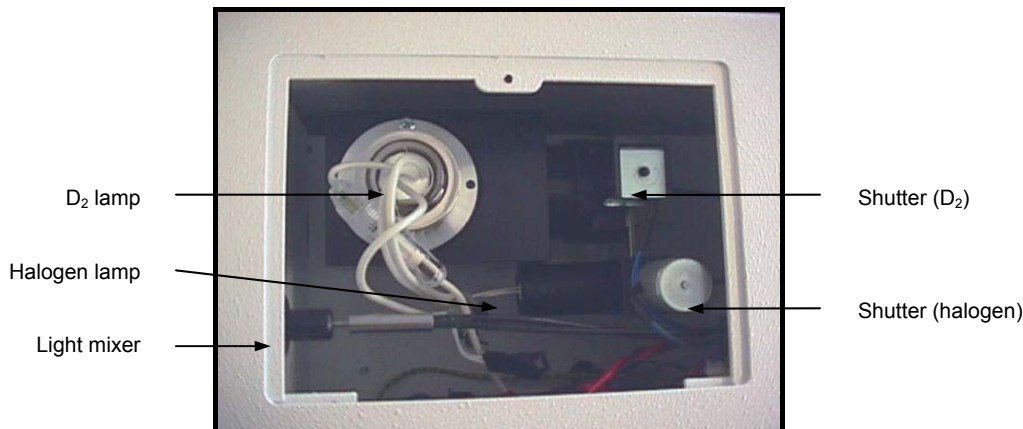
Fig. 4 Right side view (lamp replacement lid open)

The lamps are easily accessible by opening the lamp replacement lid. Also accessible are the halogen lamp, the shutter, some of the fiber optical cables and the light mixer.

THE KNAUER DAD K-2800 UV has no halogen lamp, just one shutter and no light mixer

The D₂ lamp can be removed by unscrewing the two screws, then pulling the lamp out of the lamp holder and unplugging the lamp cable. The halogen lamp is removed in the same way (just one screw). For reinstallation the lamp is put into the lamp holder, fixed with the screw(s) and afterwards the lamp cable is reconnected. The lamps are prealigned and need no further alignment.

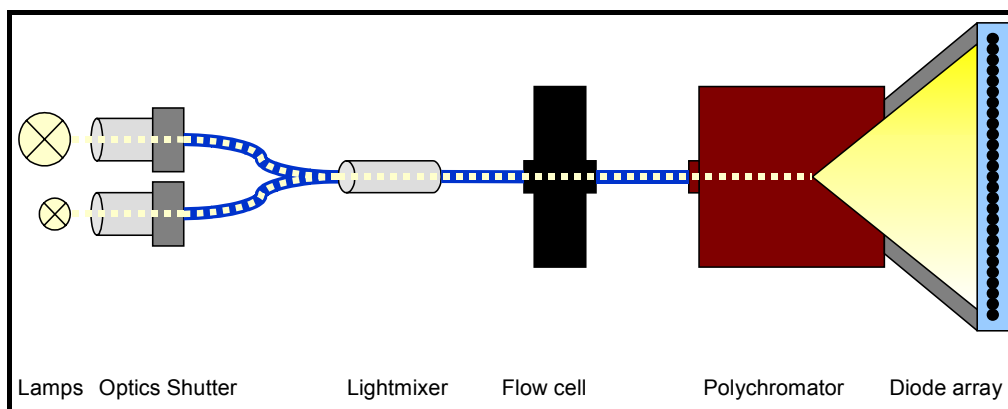
Fig. 5 *Optical system overview*

The main parts of the KNAUER Diode Array Detector K-2800 are shown in Fig. 5. The light from both lamps is focused by separate optics, passes separate shutters and goes through two fiber optical cables into the light mixer which provides perfectly balanced mixing of ultraviolet, visible and near infrared light. All light then goes through the flow cell and into the polychromator where it is split up into a spectrum of 190 to 1020 nm. The whole spectrum is detected by a 1024 diodes diode array (190-1024 nm, 0.8 nm per diode).

The KNAUER DAD K-2800 UV has no halogen lamp, just one shutter and no light mixer.

How to install DAD K-2800

PC Specifications

- 350 MHz Pentium or higher.
- 128 MB RAM (256 MB recommended) or more.
- 10 GB hard disk (20 GB recommended) or more.
- One or more PCI Slots available (PCI 2.2).
- Operating System: Win NT 4.0 SP3 or higher, Win 2000 or Win 95/98 (Win 95/98 is not recommended)
- Chromatography Software: KNAUER ChromGate system software V 2.8 or higher, KNAUER EuroChrom HPLC software Preparative V 3.01.02 or higher.

Make sure that:

- there is complete access to all devices and ports on the PC (e.g. local administrator or administrator).

- no drivers from previous DADs (e.g. K-2700) are installed.
- no energy saving mode or screensaver is active
- no virus protection system is active

Installation of the DAD K-2800

General Installation

- PC is off, power cable is not connected.
- Open PC.
- Install PCI card (supplied with DAD K-2800).
- Close PC.
- Connect power cable.
- Connect DAD K-2800 to PCI card via the RS-422 cable (supplied with DAD K-2800).
- Switch on PC.
- Enter BIOS (e.g. by pressing 'F10' key on Compaq PCs).
- Select *PCI BUS MASTER 'enabled'* if possible.
- Assign an **unshared** interrupt for the PCI card if possible (In general, the default interrupt is shared between two or more devices).
- Save settings and exit BIOS.
- Exit any automatic driver installation.

Installation for KNAUER ChromGate

- Install ChromGate, Hotfixes and ChromGate licenses.
(if ChromGate and/or licenses are already installed this step can be skipped)
- Install DAD K-2800 drivers from the ChromGate CD
(\Drivers\K2800\Setup.exe).
- Only if **Win 2000** is the operating system, the driver for K-2800 must be updated (go to the Device Manager and then click on BBK-PCI light → properties → update driver).
The correct driver for Windows 2000 can be found on the KNAUER ChromGate CD (\Drivers \K2800\Win2k\)
- Restart the PC.
- Check whether the yellow LED on the PCI card is on. If it is not on, the PCI card and/or driver have not been properly installed.
- If **Win 95/98** is the operating system, go to the Device Manager where the PCI card is defined as 'Linkadapter', 'J&M PCI Transputer Linkadapter'. Double click on 'J&M PCI Transputer Linkadapter' and then click on 'resources' to see whether there are any problems.
- If **Win NT 4.0** is the operating system, click on the Control Panel and check the devices. The PCI card should be defined as 'bbkpcil'. In the Event Viewer (menu sequence: Programs → Administrative Tools → Event Viewer) problems may be displayed.
- If **Win 2000** is the operating system, check the Device Manager where the PCI card is defined as 'OSLink', 'BBK-PCI light'. Double click on 'BBK-PCI light' and then click on resources to see whether there are any problems.

Do not install any 'older' software (e.g. for WIN 3.1) after ChromGate and drivers have been installed otherwise the system might be not stable.

Installation for KNAUER EuroChrom, Preparative Version:

- Install KNAUER EuroChrom, Hotfixes and license.
(if EuroChrom and license are already installed this step can be skipped)
- Install drivers for DAD K-2800 from the KNAUER EuroChrom CD.
(\Drivers\K2800\Setup.exe)
- Only if the operating system is Windows 2000 the driver has to be updated. (Device Manager → OSLink → bbkPCILight → properties → driver). The correct driver can be found on the EuroChrom CD
(\Drivers\K2800\Win2k\pci).
- Restart the PC.
- Check whether the yellow LED on the PCI card is on. If it is not on, the PCI card and/or driver have not been properly installed.
- If **Win 95/98** is the operating system, go to the Device Manager where the PCI card is defined as 'Linkadapter', 'J&M PCI Transputer Linkadapter'. Double click on 'J&M PCI Transputer Linkadapter' and then click on 'resources' to see whether there are any problems.
- If **Win NT 4.0** is the operating system, click on the Control Panel and check the devices. The PCI card should be defined as 'bbkpci'. In the Event Viewer (menu sequence: Programs → Administrative Tools → Event Viewer) problems may be displayed.
- If **Win 2000** is the operating system, check the Device Manager where the PCI card is defined as 'OSLink', 'BBK-PCI light'. Double click on 'BBK-PCI light' and then click on resources to see whether there are any problems.

DAD K-2800

- Switch on the DAD K-2800 (after about two seconds the red Error LED will flash once).
- If **ChromGate** software is used, the detector's lamp/lamps will be switched on when a method is opened (busy LED is on, UV LED is on after about 20 seconds, VIS LED is on if halogen lamp is installed).
- If **EuroChrom Preparative Version** software is used, the detector's lamp/s will be switched on after clicking on 'Lamp Ctrl' or 'Info' and 'Reset' in the detector's hardware setup.

Troubleshooting

If problems occur:

- Disable the sound and/or USB in the PC's BIOS.
- Deinstall the driver for the PCI-card and reinstall it from the Software CD
- Plug the PCI card into another PCI slot.

In addition, please refer to the ChromGate or EuroChrom Preparative Version manuals.

Operating the Detector with ChromGate

Detector setup

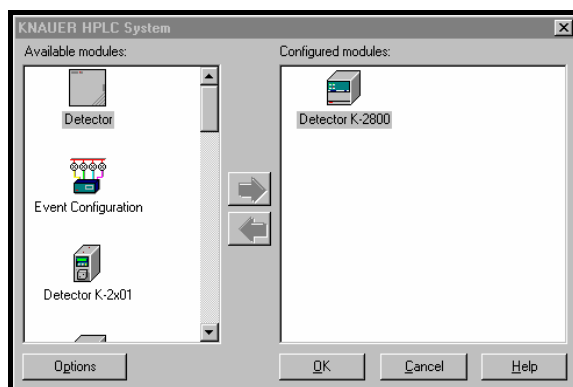
Fig. 6 KNAUER HPLC System

In the KNAUER HPLC System setup Detector K-2800 is chosen and by double clicking on the detector the configuration menu is entered.

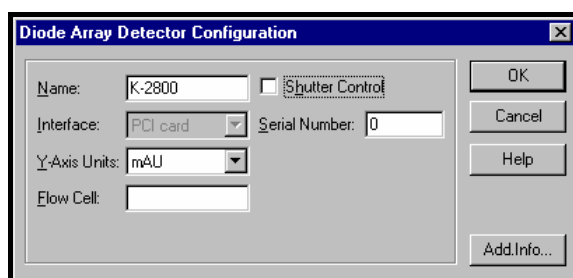
Fig. 7 Diode Array Detector Configuration

In the Diode Array Detector Configuration menu it is possible to change settings for:

- Name (default: K-2800)
- Y-Axis Units: AU, mAU or μ AU (default: mAU)
- Flow Cell: The type of installed flow cell can be entered (default: blank)
- Shutter Control: if this box is marked shutter control is enabled (default: disabled)
- Serial Number: The serial number of the detector can be entered (default: 0)

By clicking on OK the changes are saved and the Diode Array Detector Configuration menu is closed.

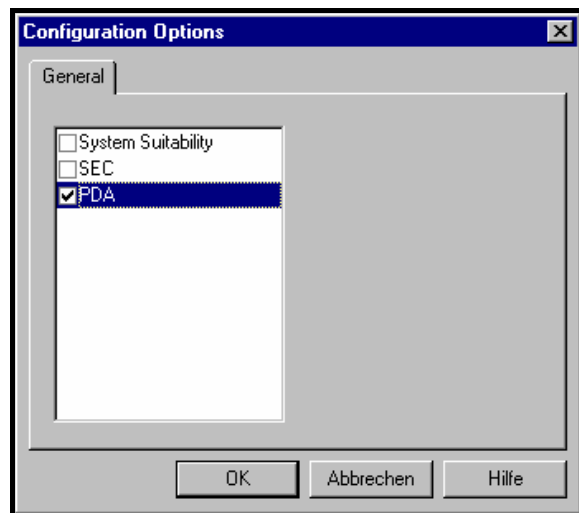
Fig. 8 Configuration Options

If one wishes to work in PDA mode (spectra acquisition enabled) the PDA option must be activated from the Options menu.

(PDA option can only be activated if a PDA license has previously been installed !)

Settings and initial tests

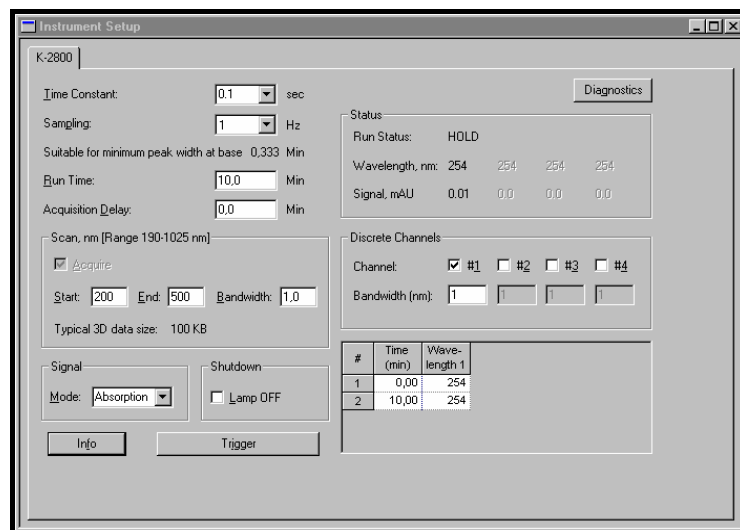
Fig. 9 Instrument Setup

When the Instrument is opened by double clicking on it's icon, ChromGate initializes the DAD K-2800 and switches on the lamp(s). (D₂-lamp needs about 20 seconds warm up time)

In the Instrument Setup menu (accessible via Method / Instrument Setup) the method's main settings can be selected.

- Time constant: 0.1, 0.2, 0.5, 1, 2, 5, 10 sec (default: 0.1 sec)
- Sampling: 0.5, 1, 2, 5 Hz (default: 1 Hz)
- Run Time: (default: 10 Min)
- Acquisition Delay: (default: 0 min)
- Scan (If the PDA option is activated): Start, End and Bandwidth can be chosen within the shown range (e.g. 190-1025 nm – default: 200-500 nm)
- Bandwidth: (default: 1 nm, recommended: 1 nm)

- Signal Mode: Absorption / Intensity (default: absorption)
- Shutdown Lamp off: if lamp off is enabled the lamp(s) will be switched off at the end of a run (default: disabled)
- Channel: up to four separate channels can be activated for data acquisition (default: one channel)
- Bandwidth: for each channel bandwidth can be set separately (default: 1 nm, recommended: 4 nm)
- Detector (Time) Program: for every channel the wavelength can be programmed separately – the entered wavelengths must be within the scan range

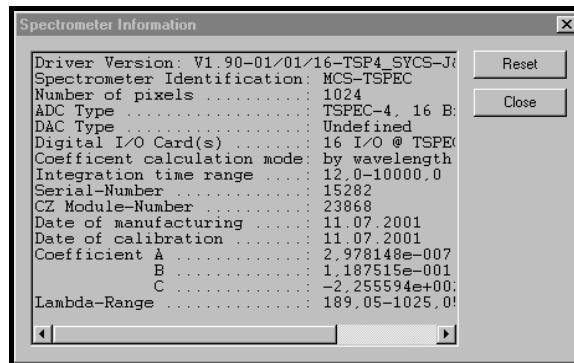
Fig. 10 Spectrometer Information

By clicking on Info and then on Reset, the Spectrometer Information from the current detector will be loaded. The individual wavelength range of this detector will now also be shown as Range in the Instrument Setup (there is no information about what lamps are installed in the DAD). All parameters are saved together with the current method. These only need to be updated if a new method is created or if the detector is exchanged.

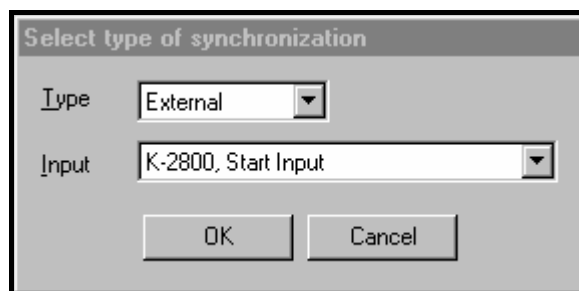
Fig. 11 Select type of synchronization

In the Trigger menu the trigger can be chosen between None (immediate start), Manual (start after confirming with Enter), External (Trigger must be applied to a chosen instrument) or Wired (Trigger must be applied to all instruments). (default: None)

Input defines the instrument which should be connected to the trigger.

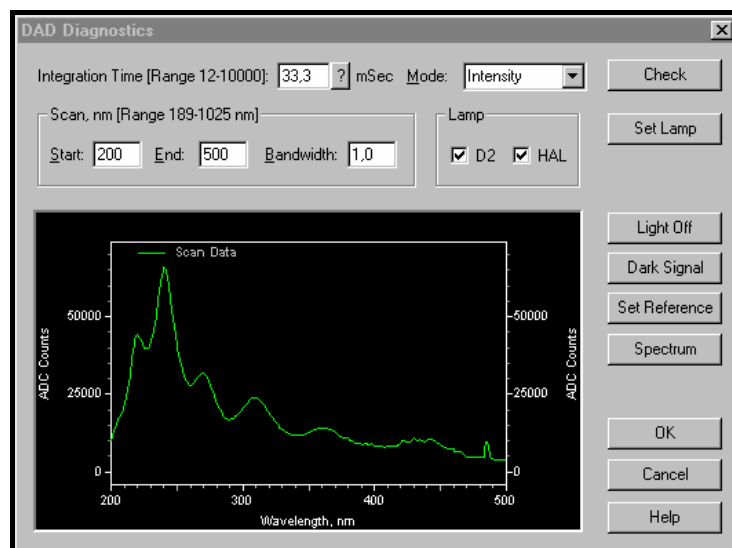
Fig. 12 DAD Diagnostics

Clicking on Diagnostics opens the DAD Diagnostics window

- Scan Range and Bandwidth are set according to the method's scan range but can be changed for diagnostics
- Integration Time: by clicking on ? the current light intensity on the diode array is measured and an optimized integration time is calculated and displayed after a few seconds.
(Settings: flow cell installed; scan range: 200-500 nm; flow rate: 1 ml/min methanol)
The Integration Time should be below 50 mSec for a completely new detector and below 100 mSec in any other case.
- Mode : Absorption / Intensity (default: Intensity)
- Lamp: on / off can be selected for D₂ and halogen lamp – afterwards Set Lamp has to be clicked on ! (default : on)

There is no information that is sent from the detector to the PC regarding whether a lamp is on or off. The PC has no information about what lamp is installed, meaning that the halogen lamp can be switched on but if it is not installed nothing happens.

- Check: This function is not yet available
- Light Off: if shutter control is activated the shutter can be opened / closed (default: closed)
- Dark Signal: Dark signal can be measured
- Set Reference: a reference spectrum can will be measured
- Spectrum: click on this button to measure an absorption or intensity

Operating the Detector with EuroChrom, Preparative Version

Detector setup

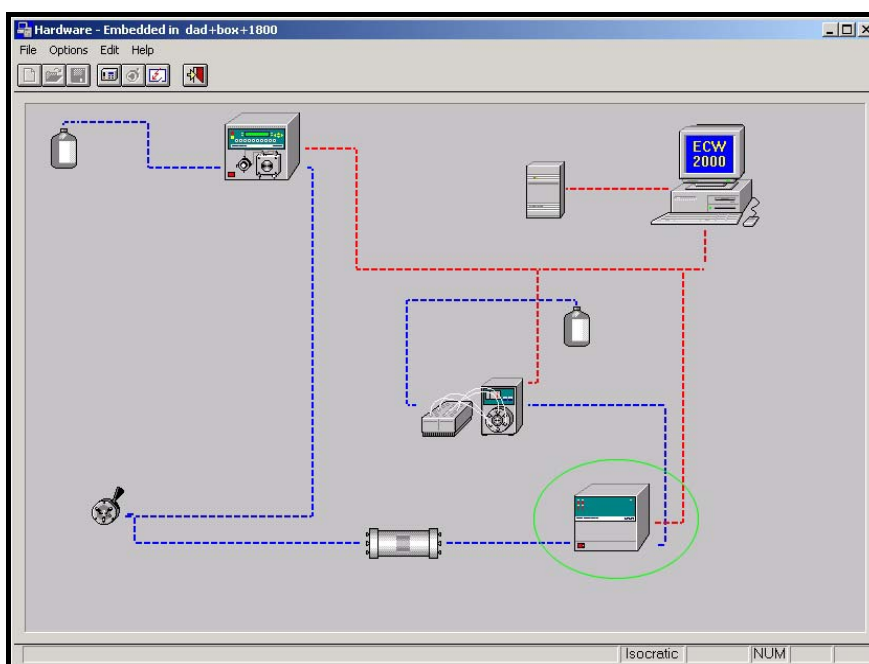
Fig. 13 Hardware

From the Hardware window, the Detector Setup menu is opened by selecting and double-clicking on the K-2800 detector icon (circled green).

The 'Detector Setup' dialog box contains the following settings:

- Name: Det1
- Channels: A B C D
- Process: A B C D
- Integration time [ms]: 6
- Scans to average: 1
- Averaging: On Off
- Wavelength range: 188 - 1023 nm
- Begin scan at: 200 nm
- End scan at: 600 nm
- Range to average: 0.8 nm
- Data: Max [mAU]:
 - Absorbance 2000
 - Emission 100
 - Transmission 100
 - ADC Counts 15000
- An. Output Scale: 10 V 5 AU (used for Absorbance only)
- Interface board type: PCI
- Shutter Control
- Comments: (empty text area)
- Buttons: OK, Cancel, Default, Lamp Cntl, Info
- Footer: Measurement with current settings will produce 0.1 MB of data.

Fig. 14 Detector Setup

In the Detector Setup the following settings can be entered:

- Name (default: Det1)
- Channels: Up to four separate channels can be selected for data acquisition (default: one channel)
- Process: To decide whether or not the measured detector signal shall be evaluated.
- Averaging: enables/disables averaging (default: enabled)
- Integration time (ms): As a rule integration time should not exceed 100 ms. Otherwise there may be a problem with the light intensity.
- Begin / End scan at: Can be chosen within the detector's individual scan range (default: 200 nm – 600 nm)
- Analog Output Scale: It can be chosen which Absorption (in AU) should equal 10 V on the analog output(s).
- Data: It can be selected whether Absorption, Emission, Transmission or ADC Counts should be recorded. (default: Absorption)
- Shutter Control: Shutter(s) can be enabled or disabled. (default: disabled)

By clicking on OK all settings are saved and the Detector Setup menu is closed.

| | |
|---------------------------|----------------|
| Owner: | KNAUER |
| Serial number: | 25292 |
| Module number: | 27048 |
| Date of manufacture: | 3 / 13 / 2002 |
| Date of calibration: | 3 / 13 / 2002 |
| Calibration coefficients: | a = +3.521e-07 |
| | b = +1.182e-01 |
| | c = -2.233e+02 |
| Integration range [ms]: | 12 - 10000 |
| Wavelength range [nm]: | 187.9 - 1023.6 |
| Pixels in array: | 1024 |

Fig. 15 Diode Array Detector

By clicking on Info the Diode Array Detector information will be shown and updated after clicking on Reset. This information is stored in the Hardware configuration after clicking on OK. Afterwards, the exact wavelength range of the current detector is shown as Wavelength range in the Detector Setup (there is no information about the installed lamp(s)). All parameters are stored with the current method and they only need to be updated for a new method or if the detector is exchanged.

Settings and initial tests

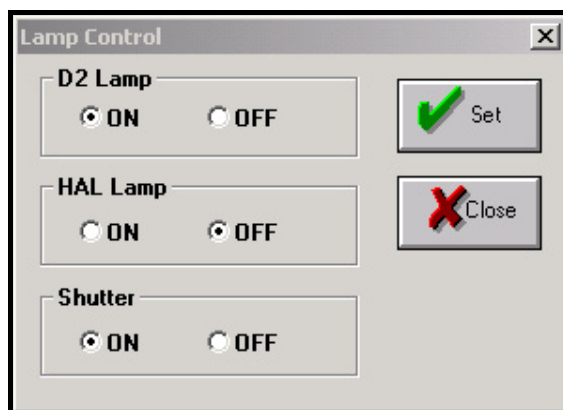
Fig. 16 Lamp Control

Clicking on Lamp Cntl in the Detector Setup opens the Lamp Control menu where the lamps and shutter(s) can be switched on and off.

Lamp: on / off can be set separately for D₂ and halogen lamps – afterwards Set must be clicked to switch the lamps and shutter asselected ! (default: on)

Only from the detector's operation display, can the user see which lamp is on and which lamp is off. The software receives no information about the lamp status.

PC and software can not 'see' what lamp(s) is (are) installed in the DAD. For example a halogen lamp can be switched on and off through the software although it might not actually be installed.

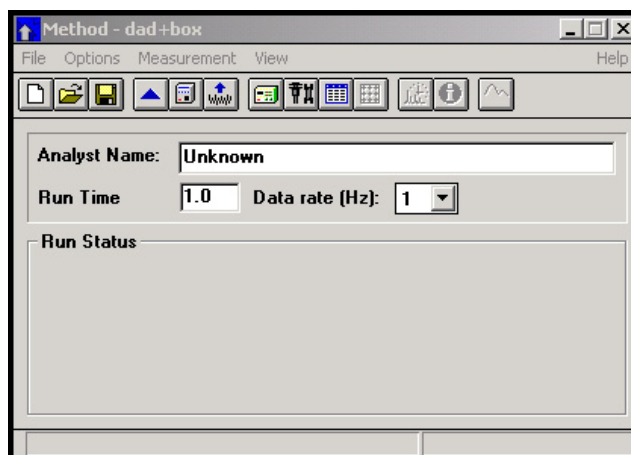
Fig. 17 Method

From the Method menu the user can access all other important menus. Also Run Time and Data rate (Hz) can also be set here.

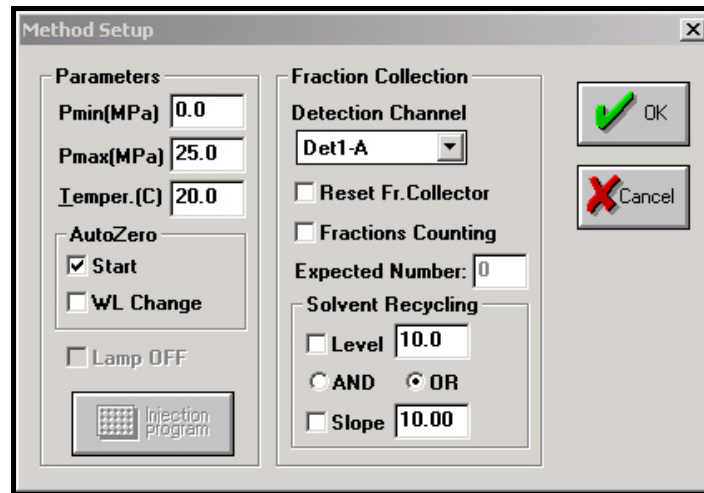
Fig. 18 Method Setup

In the Method Setup menu Auto Zero and the Detection Channel for Fraction Collection can be configured.

| | Time min | Flow ml/min | Pump %A | D.In | WLength 1-A |
|---|----------|-------------|---------|------|-------------|
| 1 | 0.00 | 1.000 | 100 | - | 254 |
| 2 | 1.00 | 1.000 | 100 | - | 254 |

Fig. 19 Time table

From the Time table menu a wavelength program can be set for every enabled channel.

Spare Parts and Accessories:

| Article | Order No |
|---|----------|
| Upgrade + Upgrade Kit from UV instrument to UV/VIS/NIR instrument | A4446 |
| deuterium spare lamp | A4447 |
| halogen spare lamp | A4448 |
| Analytical flow cell, 10 mm, VA, 1/16", 20 ml/min, 300 bar | A4130 |
| Analytical flow cell, 3 mm, VA, 1/16", 50 ml/min, 300 bar | A4131 |
| Analytical flow cell, 3 mm, PEEK, 1/16", 10 ml/min, 30 bar | A4132 |
| Preparative flow cell, 0.5/1.25/2 mm, VA, 1/8", 1,000 ml/min, 200 bar | A4133 |
| Preparative flow cell, 0.5/1.25/2 mm, PEEK, 1/8", 1,000 ml/min, 100 bar | A4134 |
| Preparative flow cell, 0.5/1.25/2 mm, VA, 1/4", 10,000 ml/min, 200 bar | A4135 |
| Preparative flow cell, 0.5 mm, VA, 1/16", 250 ml/min, 200 bar | A4136 |
| Preparative flow cell, 0.5 mm, PEEK, 1/16", 250 ml/min, 100 bar | A4137 |
| Holmium oxide test cell | A4129 |

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Hegauer Weg 38
14163 Berlin, Deutschland

**WellChrom Diode Array Detector K-2800,
Order number A 4444, A 4445**

complies with the following requirements and product specifications:

- Low Voltage Ordinance (73/23/EWG);
EN 61010 – 1 (1993)
- Engineering Guidelines (89/392/EWG)
- EMV Ordinance (89/336/EWG)
EN 50081 – 1 (1992)
EN 55011 (1991) Class B
EN 55022 (1987) Class B
EN 50082 – 1 (1992)
IEC 801 – 2 (1984),
includes IEC 41 B (sec) 81 (1992)
IEC 801 – 3 (1984)
IEC 801 – 4 (1988)

The product was tested in a typical configuration.

Berlin December, 10th 2002
Bernward Rittgerodt (Managing Director)

The CE Shield is attached to the rear of the instrument.

Guarantee statement

The guarantee period of the WellChrom Diode Array Detector K-2800 is 12 months beginning from the date of dispatch from Berlin. Operation inconsistent with manufacturer's instructions or damage caused by unauthorized service personnel are excluded from guarantee. Damage caused by blockages and wear and tear parts such as fuses, lamps, and fiber optical cables are not covered by the guarantee. Claims under this guarantee are valid only if the enclosed guarantee card is returned to us at the address shown below within 14 days of receipt of the instrument. Defective detectors should be sent to the manufacturer for repair.

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If we find a defect covered by the guarantee, repair or replacement, at our discretion, will be carried out free of charge. Packing and transport costs are borne by the purchaser.

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Dr. Ing. Herbert Knauer GmbH