
Instructions for Adjusting the Path Length in Preparative Flow Cells

V4060 06/2001

These instructions are valid for the preparative flow cells A4066, A4067, and A4068 with 1/8" or 1/4" connectors.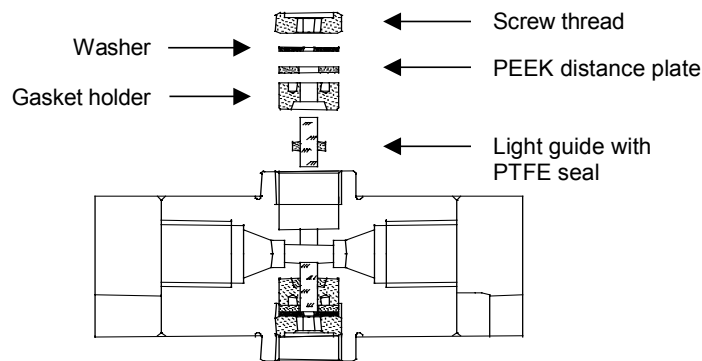

Sectional view of a preparative flow cell

The measurement path length is set to 2mm when manufactured. The measurement path length can be adjusted to 2; 1.25 or 0.5 mm. If desired, follow the below steps to reduce the path length to 1.25 or 0.5 mm

1. Loosen the cap nut with a hexagonal head wrench.
2. Remove the stainless steel sealing ring and the PEEK spacer.
3. Replace the stainless steel ring and retighten the nut.
Without a spacer the light flow will be pushed deeper into the flow cell. (0,75 mm), resulting in a reduction in the measurement path length. To reduce the measurement path length even further to 0,5 mm use the other side of the cell in analog mode.
- To increase the measurement path length back to 0,75 mm reinsert the spacer.
4. Loosen the cap nut, remove the stainless steel sealing ring, and remove the light guide mounting with a pair of tweezers.
5. Use a clean towel to push the light guide, about 1 mm, outwards in order to increase the path length. Avoid touching the light flow with fingers.
6. Reinsert the mounting in the cell.
7. Replace the PEEK spacer and then the stainless steel ring.
8. Carefully retighten the cap nut.

Tightening the cap nut will push the rod shaped light guide into the right position. Inserting the spacer increases the path length to 0.75 mm. When changing the path length it is not necessary to change the PTFE seal.

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